

# EnviroMon Data Generator




The EnviroMon Data Generator allows you to extract data by a few simple clicks and without having to go into the software.

In order for the software to work, you must place it in the same location as your EnviroMon folder. By default this can be found in:

C:\Program Files\Pico Technology\EnviroMon\

You may also need to download the .NET 2 framework, as the program is written in C#. The program works with Windows XP and Vista.

The image below shows the buttons available and what they all mean:

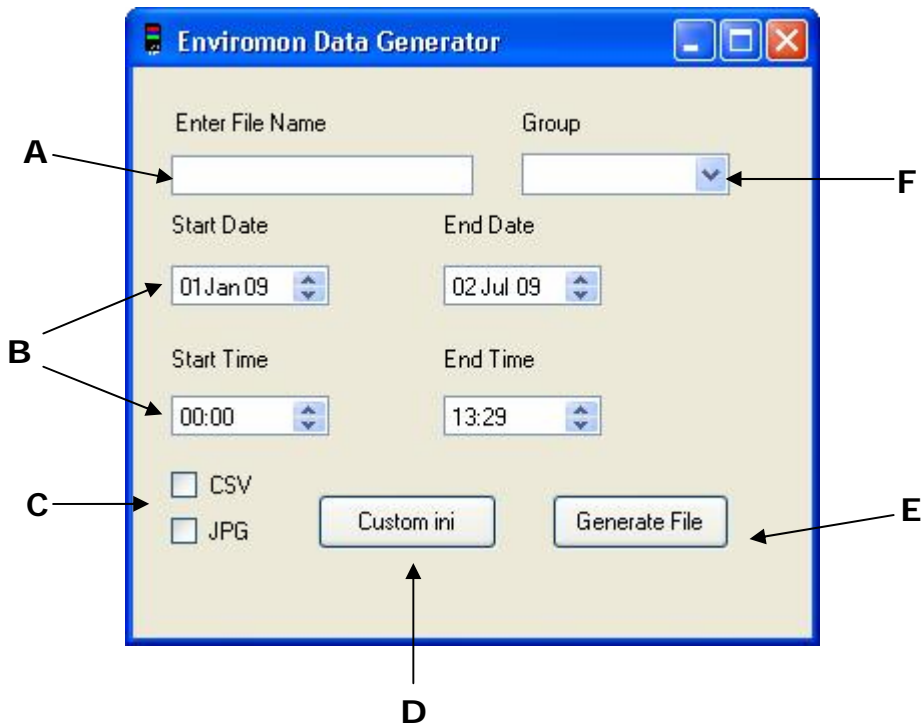

- A. Enter the name of a file you want to create.
- B. Enter time range of data you want to view.
- C. Choose the format of the file, either a .csv, .jpg or both.
- D. If you have a .ini file located in a different location, browse to its location.
- E. Once all options are chosen, "Generate File" creates the file(s).
- F. If you have created sensor groups in EnviroMon, all the group names will appear in this list. Choose the group that you want to view.

[www.picotech.com](http://www.picotech.com)